

# INSTRUCTIONS BX43 SYSTEM MICROSCOPE

This instruction manual is for the Olympus System Microscope Model BX43. To ensure the safety, obtain optimum performance and to familiarize yourself fully with the use of this microscope, we recommend that you study this manual thoroughly before operating the microscope. Retain this instruction manual in an easily accessible place near the work desk for future reference.





In accordance with European Directive 2002/96/EC on Waste Electrical and Electronic Equipment, this symbol indicates that the product must not be disposed of as unsorted municipal waste, but should be collected separately.

Refer to your local Olympus distributor in EU for return and/or collection systems available in your country.

**NOTE:** This equipment has been tested and found to comply with the limits for a Class A digital device, pursuant to Part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his own expense.

**FCC WARNING:** Changes or modifications not expressly approved by the party responsible for compliance could void the user's authority to operate the equipment.

# **CONTENTS**

Correct assembly and adjustments are critical for the microscope to exhibit its full performance. If you are going to assemble the microscope yourself, please read section 10, "ASSEMBLY" (pages 33 to 37) carefully.

IMP	<b>ORTANT</b> – Be sure to read this section for safe use of the equipment. – 1-3
1	MODULE NOMENCLATURE 4
2	CONTROLS 5-7
3	FLOW OF OBSERVATION 8,9
4	SIMPLIFIED OBSERVATION PROCEDURE 10-12
	4-1 Basic Operation (Until Observation of Specimen)
	4-2 Microscope Adjustments (How to improve the Observed image)
	Adjusting the Interpupillary2 Adjusting the Left and RightDistance
	3 Adjusting the Centering12 4 Adjusting the Contrast
5	USING THE CONTROLS 13-25
	5-1 Base
	Adjusting the Brightness
	3 Using the Filters (Halogen Bulb Operation Only)
	5-2 Focusing Block 16,17
	Focusing Controls 16 2 Adjusting the Focus 16
	MODULE NOMENCLATURE         CONTROLS       5         FLOW OF OBSERVATION       8         SIMPLIFIED OBSERVATION PROCEDURE       10-         4-1 Basic Operation (Until Observation of Specimen)       10;         4-2 Microscope Adjustments (How to improve the Observed image)       11;         I Adjusting the Interpupillary       I Adjusting the Left and Right         Distance       11         I Adjusting the Centering       12         I Adjusting the ContrROLS       13-2         S-1 Base       13-1         I Adjusting the Brightness       13         I Adjusting the Brightness       13         S Using the Filters (Halogen Bulb Operation Only)       14         5-2 Focusing Block       16         I Adjusting the Fine Adjustment Knob       17         G Replacing the Fine Adjustment Knob       17         G Pre-focusing Lever       17         5-3 Stage       18         I Placing the Specimen       18         I Adjusting the X- and Y-Axis Knob Tension       19         G Adjusting the Stage Height       20         5-4 Observation Tube       21         I Adjusting the Diopter       21         I Adjusting the Eypejece Micrometer Disk       21
	с. С
	5 Adjusting the Stage Height 20
	5-4 Observation Tube
	Adjusting the Diopter
	5 Adjusting the Tilt
	5-5 Condenser

	5-6 Immersion Objectives	
	1 Using Immersion Objectives 25	
	5-7 Objectives with Correction Collar	
6	CAMERA RECORDING	26
	System Chart 26 2 Selecting the Camera Adap Magnification	
7	TROUBLESHOOTING GUIDE	27-29
8	SPECIFICATIONS	30,31
9	<b>OPTICAL CHARACTERISTICS</b> (UIS2 Series)	32
10	ASSEMBLY – See this section for the replacement of the light bulb. –	33-37
11	HALOGEN LAMP SOCKET INSPECTION SHEET	38
	PROPER SELECTION OF THE POWER SUPPLY CORD	39,40

# **IMPORTANT**

# A SAFETY PRECAUTIONS

- (Fig. 1)
- 1. After the equipment has been used in an observation of a specimen that is accompanied with a potential of infection, clean the parts coming in contact with the specimen to prevent infection.
  - Moving this product is accompanied with the risk of dropping the specimen. Be sure to remove the specimen before moving this product.
  - In case the specimen is damaged by erroneous operation, promptly take the infection prevention measures.
  - Follow the procedures described in Chapter "Getting Ready" (see on page 2) prior to using the accessories of this product. Otherwise, the stability of the device will be lost and the dropped specimen will cause the possibility of infection.
- 2. Install the microscope on a sturdy, level table or bench.
- 3. Be careful not to tilt the microscope too much. Otherwise, the mobile parts such as the light path selector knob may move in unintended directions.
- 4. Always use the power cord provided by Olympus. If no power cord is provided, please select the proper power cord by referring to the section "PROPER SELECTION OF THE POWER SUPPLY CORD" at the end of this instruction manual.

If the proper power cord is not used, product safety performance cannot be warranted.


5. Always ensure that the **grounding terminal** of the microscope and that of the wall outlet are properly connected. If the equipment is not grounded, Olympus can no longer warrant the electrical safety performance of the equipment.

## Operation Using the LED Lamp

Safety Note on LED (Light Emitting Diode)

The LED incorporated in the LED lamp housing is a class 1 LED product as defined in the applicable standard. The light emitted by LED is basically safe, but do not look at the illumination light <u>directly for an</u> extended period to prevent your eye from being injured

## CLASS 1 LED PRODUCT

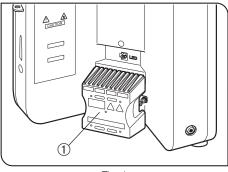

#### Fig. 1

## Operation Using the Halogen Bulb

1. To avoid potential shock hazards and burns when replacing the halogen bulb, set the main switch of the TL4 power supply unit to " **O**" (OFF) and allow the lamp socket ① and bulb to cool before touching them.

Designated bulb

## 6V30WHAL (PHILIPS 5761)

## CAUTION

# Always use the designated lamp bulb. Using other bulb may cause malfunction.

- 2. If the lamp cable or power cord comes in contact with the lamp socket or its surroundings, the cable or cord may be melt by the heat, causing an electric shock or fire hazard. To prevent this, distribute the lamp cable and power cord at enough distances from the lamp socket.
- 3. The standard service life of the lamp housing is <u>eight (8) years</u> of use <u>or</u> <u>20,000 hours</u> of total power ON period, <u>whichever is the shorter period</u>. For details, see Inspection Sheet on page 38.

## Safety Symbols

The following symbols are found on the microscope. Study the meaning of the symbols and always use the equipment in the safest possible manner.

Symbol	Explanation			
	Indicates that the surface becomes hot, and should not be touched with bare hands.			
	dicates a non-specific general hazard. Follow the description given after this symbol or in struction manual.			
I	Indicates that the main switch is ON.			
0	Indicates that the main switch is OFF.			

## Caution engraving/label

Caution engraving and label are placed at parts where special precaution is required when handling and using the microscope. Always heed the cautions.

Caution engraving position	Halogen lamp socket [Caution against high temperature]	
Caution label position	Rear panel [Caution against high temperature]	

If a caution engraving or label is dirty or peeled off, contact Olympus for the replacement or other inquiry.

## Getting Ready

1

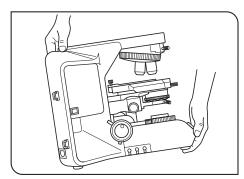


Fig. 2

- A microscope is a precision instrument. Handle it with care and avoid subjecting it to sudden or severe impact. Also note that the microscope does not have a waterproof construction.
- 2. Do not use the microscope where it is subjected to direct sunlight, high temperature and humidity, dust or vibrations. (For the operating conditions, refer to section 8, "SPECIFICATIONS".)
- 3. When moving the microscope, remove the observation and specimen and carefully carry the microscope by holding the grasping part on the rear of the arm and the front part of the base as shown in Fig. 2.

CAUTION

- To prevent damage to the microscope, never hold it by the stage or coarse/fine adjustment knob.
- Even when moving the microscope along the desktop surface, be sure to remove the specimen in advance.
- When the microscope needs to be packaged for forwarding to a remote location, contact Olympus.
- 4. When the accessories of this product are used, make sure that the system height does not exceed 1 meter. Otherwise, the stability of the device will be lost.

## 2 Maintenance and Storage

1. To clean the lenses and other glass components, simply blow dirt away using a commercially available blower and wipe gently using a piece of cleaning paper (or clean gauze).

If a lens is stained with fingerprints or oil smudges, wipe it gauze slightly moistened with commercially available absolute alcohol.

CAUTION • Since the absolute alcohol is highly flammable, it must be handled carefully.

Be sure to keep it away from open flames or potential sources of electrical sparks — for example, electrical equipment that is being switched on or off, which could cause ignition of a fire. Also remember to always use it only in a well-ventilated room.

- Set the main switch to " **O** " (OFF) and wait for the lamp housing to cool down sufficiently before proceeding to maintenance.
- 2. Do not use organic solvents to clean the microscope components other than the glass components. To clean them, use a lint-free, soft cloth slightly moistened with a diluted neutral detergent.
- 3. Do not disassemble any part of the microscope except for the parts that are specified to be disassembled in this manual, as this could result in malfunction or reduced performance.
- 4. After using the microscope, set the main switch to "**O**" (OFF), (wait for the lamp socket to cool down sufficiently if the halogen bulb has been used,) and keep it covered with a dust cover during storage.
- 5. Before disposing of this product, be sure to follow the regulations and rules of your local government.

# 3 Caution

If the microscope is used in a manner not specified by this manual, the safety of the user may be imperiled. In addition, the equipment may also be damaged. Always use the equipment as outlined in this instruction manual.

\_\_\_\_\_

The following symbols are used to set off text in this instruction manual.

CAUTION : Indicates a potentially hazardous situation which, if not avoided, may result in minor or moderate injury or damage to the equipment or other property. It may also be used to alert against unsafe practices.
 Indicates commentary (for ease of operation and maintenance).

## 4 Intended use

This instrument has been designed to be used to observe magnified images of specimens in various routine work and research applications.


Do not use this instrument for any purpose other than its intended use.

This instrument is considered a medical device in the following countries: US, and China.




This device complies with the requirements of directive 98/79/EC concerning in vitro diagnostic medical devices. CE marking means the conformity to the directive.


The modules mentioned below show only the typical product names. As there are some products that are not mentioned but also applicable to this microscope, check the latest catalogues or consult Olympus. For the products marked " \* ", also read their instruction manuals.

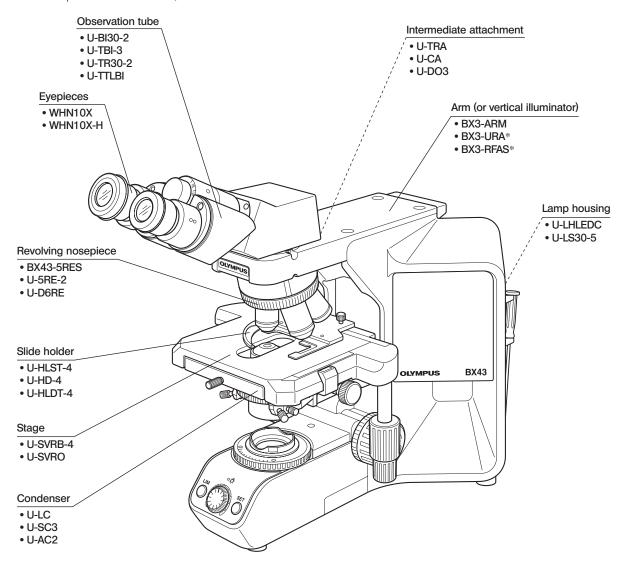





Olf you have not yet assembled the microscope, read section 10, "ASSEMBLY" (pages 33 to 37).

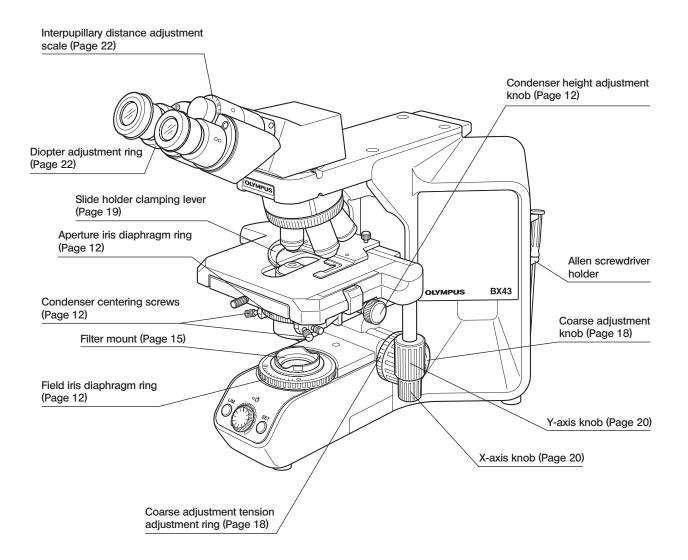

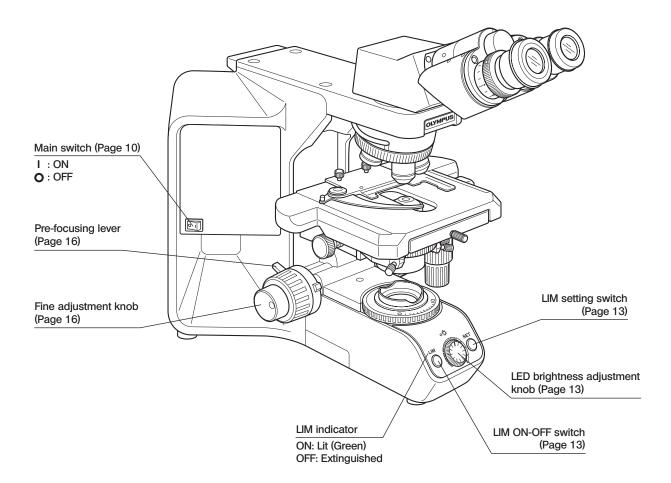

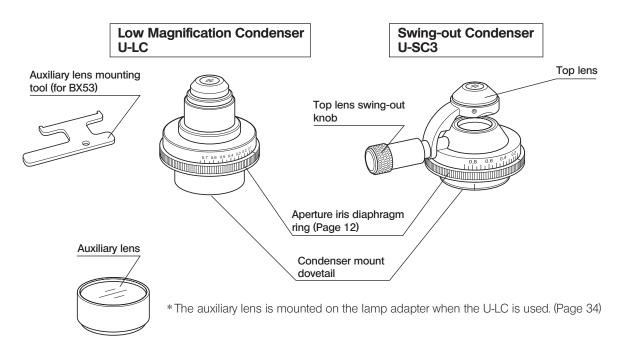


# «Modules for halogen lamp operation»

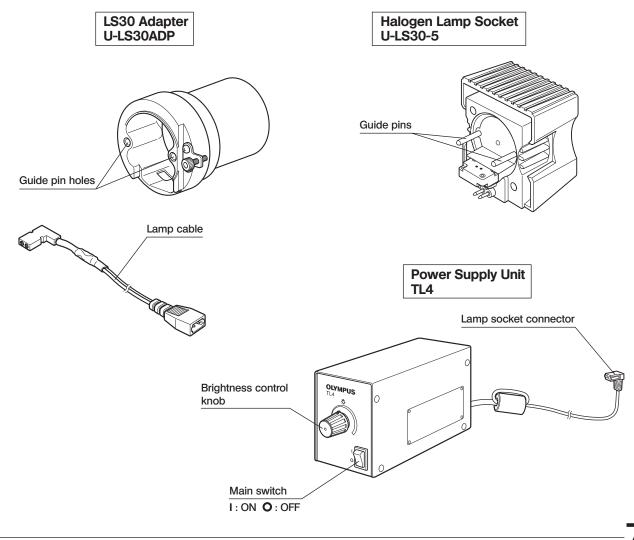


# **FLOW OF OBSERVATION**

When the LED lamp is used and the LIM is set, the LED brightness adjustment knob is defeated.When the halogen bulb is used, set the LBD filter in the filter mount on the base (or in the U-FC filter cassette).

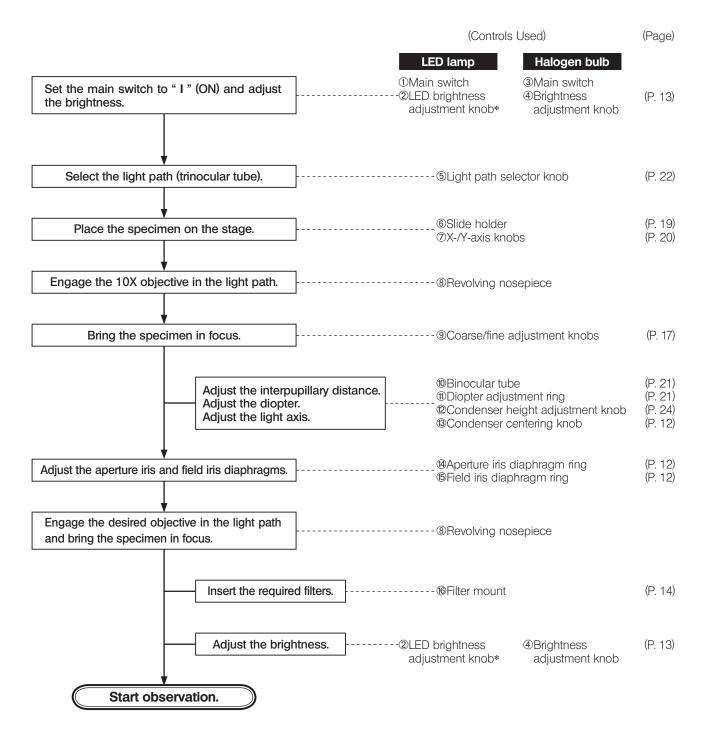

\* The LED brightness adjustment knob is defeated when the LIM is set (see page 34).

**BX43** 

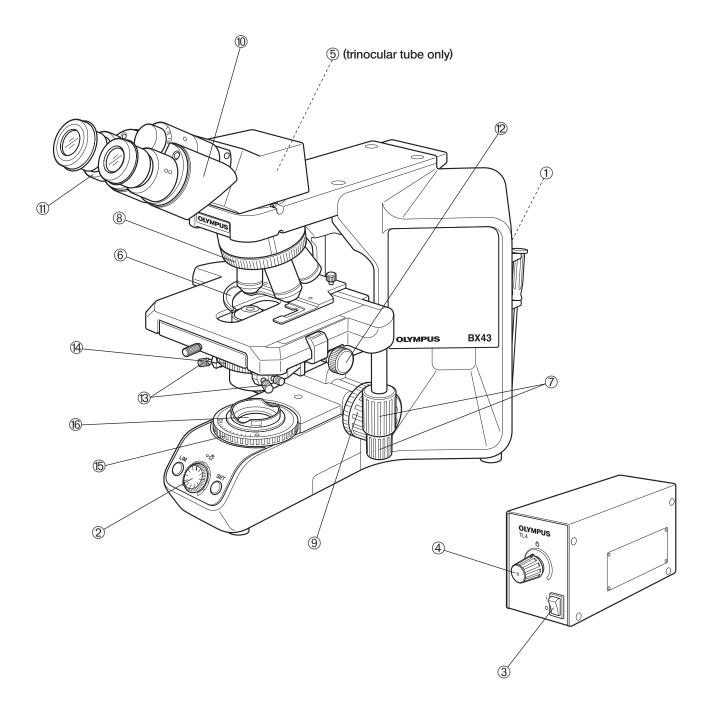



©Copy the observation procedure pages on separate sheets and post it near your microscope.

# **4** SIMPLIFIED OBSERVATION PROCEDURE

# 4-1 Basic Operation (Until Observation of Specimen)

This section describes the basic operation of the microscope until the start of observation of a specimen. For the detailed operating procedure of each control, please read the description page specified below.

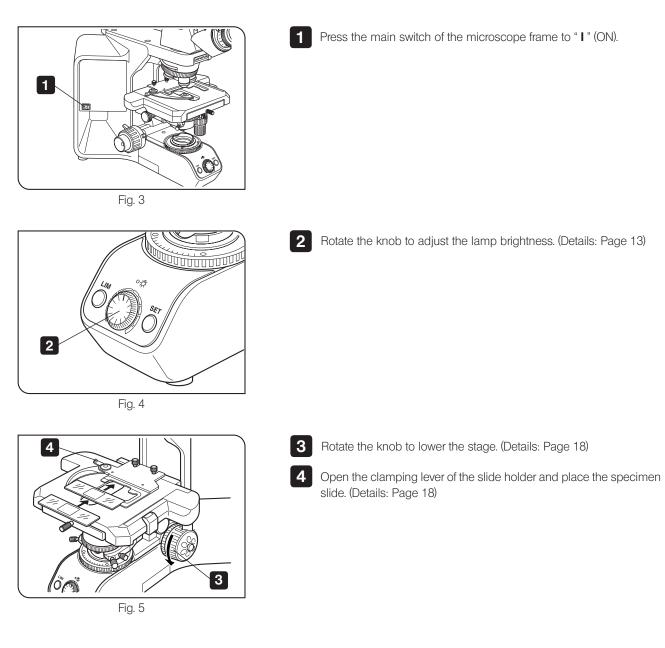

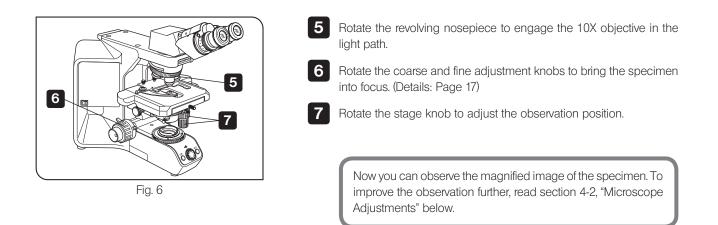


# 4-2 Microscope Adjustments (How to Improve the Observed Image)

Adjust the microscope as described below to improve the observed image.

1

2

1

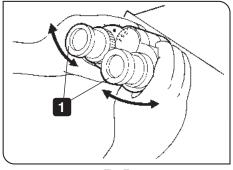

#### Fig. 7

# Adjusting the Interpupillary Distance

While looking through the eyepieces, adjust for binocular vision until the left and right fields of view coincide completely. ©Note your interpupillary distance so that it can be quickly duplicated.



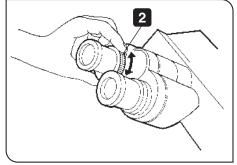


Fig. 8

# Adjusting the Left and Right Focusing (diopter adjustment)

- Looking through the eyepiece without the diopter adjustment ring, rotate the coarse and fine adjustment knobs to bring the specimen into focus.
- 2 Looking through the eyepiece with the diopter adjustment ring, turn only the diopter adjustment ring to focus on the specimen.
  - ♥If the diopter adjustment rings are provided on both the left and right sides, see "Adjusting the Diopter" on page 21.
    - The diopter adjustment rings may be provided on either the eyepieces or the observation tube.

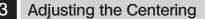

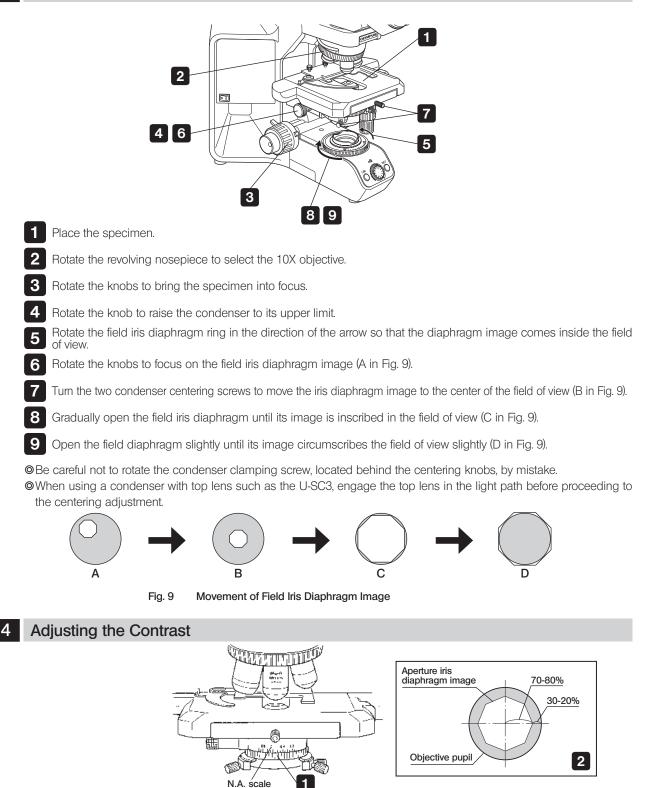


Set the condenser aperture iris diaphragm ring so that the condenser NA scale position indicates 70% of the NA of the objective in use. After this, observe the image and fine adjust the aperture iris diaphragm to the desired condition.

2 The image of the aperture iris diaphragm can be viewed by removing the eyepieces and looking into the observation tube. Adjust so that the image looks approximately like that shown in the figure above.

12

# **5** USING THE CONTROLS

# 5-1 Base

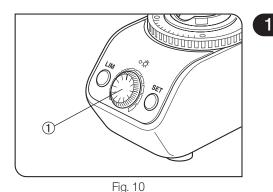

# Adjusting the Brightness

(Fig. 10)

- 1. When the LED lamp is used for observation, turn the LED brightness adjustment knob ① clockwise to make illumination brighter.
- 2. When the halogen bulb is used for observation, adjust the brightness adjustment knob (2) of the TL4 power supply unit.

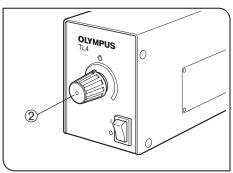


Fig. 11

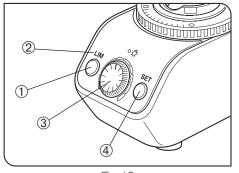


Fig. 12

## 2 Setting the LIM Brightness

(Fig. 12)

- ©LIM stands for Light Intensity Manager. It is effective when the BX43-5RES revolving nosepiece with sensor and the U-LHLEDC LED lamp housing are used.
- 1. Press the LIM ON-OFF switch 1 to "OFF" (so that the LIM indicator 2 is extinguished).
- 2. Engage a low-power objective in the light path and bring the specimen in focus.
- 3. Rotate the brightness adjustment knob ③ for optimum brightness for observation.
- 4. Press the LIM setting switch ④ to enter the current brightness.
- 5. Set the optimum brightness for the mounted objectives by repeating the above for each of them.

## To modify the LIM setting:

Engage the objective you want to set the new LIM setting in the light path, set the LIM ON-OFF switch to "OFF," and perform steps 3 and 4 above.

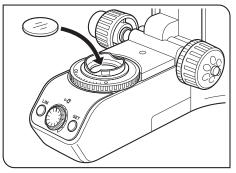


Fig. 13

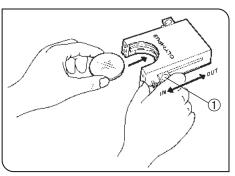


Fig. 14

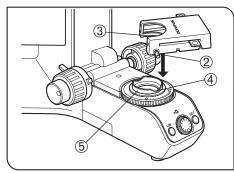


Fig. 15

# 3 Using the Filters (Halogen Bulb Operation Only)

You can place a filter in the light path with either method.

- Place a filter on the filter mount on the base and engage in the light path. (Page 14)
- Insert a filter in the U-FC filter cassette, attach it on the filter mount and engage it in the light path by sliding the filter level. (Page 14)

#### Mounting a Single Filter (Fig. 13)

A filter with a diameter of 45 mm can be placed on the filter mount on the base. If you need to place multiple filters, please purchase the U-FC filter cassette.

Even when a filter cassette is used, a filter with thickness of no more than 3 mm can still be placed on the filter mount.

#### Using the Filter Cassette (Figs. 14 to 17)

#### Loading Filters into Filter Cassette

- The filter cassette accommodates filters with a diameter of 45 mm and thickness of 2.7 mm or less.
- ©The filter cassette has two filter levels on the right side and one on the left side.
- When the filter cassette is used, the ECO mode cannot be used because the human sensor is hidden by the filter cassette.
- 1. Move all filter levers to the OUT position except for the one belonging to the slot into which the filter is to be inserted.
- 2. Slide lever  ${\rm \textcircled{O}}$  to the IN position. Make sure that it clicks securely into place. (Fig. 14)
- 3. Holding the lever in the position shown in Fig. 14, put the filter into the cassette by inserting it in the direction indicated by the arrow.
- 4. Insert the other two filters in the same manner.

#### Mounting the Filter Cassette

- 1. Fully loosen the filter cassette clamping screw 2. (Fig. 15)
- 2. Align the key ③ on the bottom surface of the filter cassette with the positioning slot ④ on the filter mount, then snap the filter cassette into place from above.
- 3. Rotate the filter cassette to align its sides with the base. (Fig. 16)
- 4. Align the clamping screw ② with the positioning hole ⑤ on the filter mount, then tighten the screw to fasten the filter cassette.



When the filter cassette is installed, the condenser may hit it when the stage is lowered. Therefore, exercise caution when lowering the stage with the filter cassette installed.

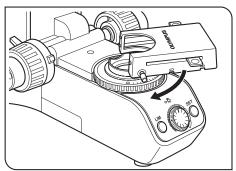


Fig. 16

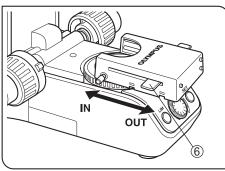


Fig. 17

## Using the Filter Cassette (Fig. 17)

Usable Filters	Applications		
45LBD-IF*	Color balancing filter		
45-ND6, 45-ND25	Neutral density filter		
45G-530, 43IF550-W45*	Green		
450-560	Orange	B&W contrast filters	

Up to three of the above filters can be loaded in the filter cassette. Moving the levers (6) on the left and right sides of the cassette to the IN position brings the corresponding filter into the light path.

- \* When an IF (interference) type filter is used, flare and ghost may be observed. These phenomena may be reduced by adding ND filter(s) to the IF filter.

## What is LBD filter? —

The color of the sun at the noon of a clear day is called the daylight color, which looks white to human eyes. On the other hand, the halogen bulb used as the light source of this microscope does not emit the daylight-color light so the LBD filter is engaged in the light path to convert the color of halogen bulb light into daylight. In general, engaging the LBD filter makes it possible to observe specimens in natural color tones.

# 5-2 Focusing Block

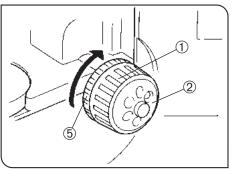


Fig. 18

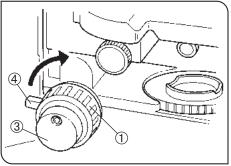



Fig. 19

# **Focusing Controls**

1

The focusing block consists of the controls listed in the following table, which also describes the functions of the controls.

Ν	lo.	Name	Function
(	1	Coarse adjustment knob	Moves the focus position widely.
(	2	Fine adjustment dial	Fine adjusts the focus position.
(	3	Fine adjustment knob	Fine adjusts the focus position. This knob can be attached to the fine adjustment dial on either the left or right side. (When the microscope is shipped, it is attached to the right-side dial.)
(	4)	Pre-focusing lever	Sets the upper limit of the stage position (to prevent contact between the speci- men and objective).
(	5	Tension adjustment ring	Adjusts the force required to rotate the coarse adjustment knob.

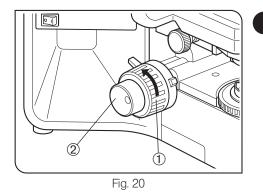

2 Adjusting the Focus

With both the coarse adjustment knob 1 and fine adjustment knob 2, rotating the knob in the direction of the arrow raises the stage (lets the specimen approach the objective).

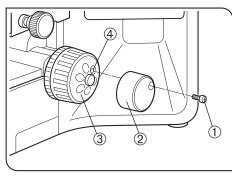



Fig. 21

# 3 Replacing the Fine Adjustment Knob (Fig. 21)

# CAUTION

# The fine adjustment knob has been attached on the right side at the factory.

- O The fine adjustment knob is designed detachable to prevent interference with hand during manipulation of the X-and Y-axis knobs.
- Usually attach the knob on the opposite side to the X- and Y-axis knobs.
- 1. Using the Allen screwdriver, loosen the clamping screw ① and remove the fine adjustment knob ②.
- 2. Remove the seal from the fine adjustment knob screw hole on the other side and attach the knob by reversing the removal procedure.
- 3. Attach a provided seal on the screw hole ④ of the removed fine adjustment knob ③.
- The fine adjustment dial ③ can be operated with your fingertip or finger surface at the same time as manipulating the X- and Y-axis knobs.

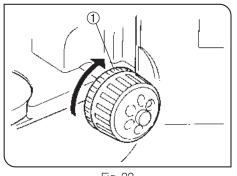

#### Fig. 22

## Adjusting the Coarse Adjustment Knob Tension (Fig. 22)

# CAUTION

Δ

# Adjust the coarse adjustment knob tension using the tension adjustment ring ①.

The coarse adjustment knob tension is preadjusted for easy use. However, if desired, you can change the tension using the tension adjustment ring (D. Turning the ring in the direction of the arrow increases tension, and vice versa.

The tension is too low if the stage drops by itself of focus is quickly lost after adjustment with the fine adjustment knob. In this case, turn the ring Fig. 22 in the direction of the arrow to increase tension.

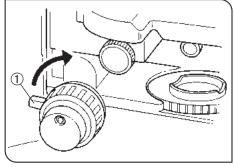


Fig. 23

## 5 Pre-focusing Lever

(Fig. 23)

The pre-focusing lever ensures that the objective does not come in contact with the specimen and simplifies focusing.

After focusing on the specimen with the coarse adjustment knob, turn this lever ① in the direction of the arrow and lock; the upper limit on coarse adjustment movement is set at the locked position.

After changing a specimen, approximate refocusing is easily accomplished by rotating the coarse adjustment knob to reach the pre-focused position. After this, all you have to do is make fine adjustments with the fine adjustment knob.

Stage movement with the fine adjustment knob is not locked.



When the pre-focusing lever is locked, a coarse movement stroke is reduced due to the microscope's mechanism, causing the stage not to be able to be lowered to the lower limit. To lower the stage to the lower limit, release the pre-focusing lever.

# 5-3 Stage

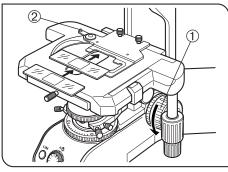


Fig. 24

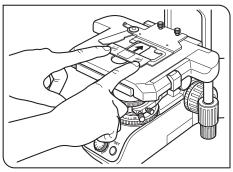


Fig. 25

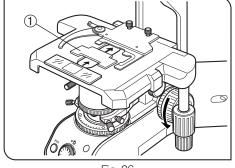


Fig. 26

# Placing the Specimen

1

2

© The dimensions of the slide glass should be 26 x 76 mm with thickness of 0.9 to 1.2 mm, and the cover glass should have thickness of 0.17 mm.

When observing very large specimens, remove the slide holder and place the specimen directly on the stage.

## Microscopy with Double-Slide Holder (Fig. 24)

1. Turn the coarse adjustment knob ① to lower the stage.

- 2. Open the spring-loaded curved finger ② on the slide holder and place one or two specimen slides on the stage from the front.
- 3. After placing the sides as far as they will go, gently release the curved finger.

## Microscopy with Single-Slide Holder (Fig. 25)

The specimen side can easily be placed by sliding it into the slide holder from the front.

## Examining a whole specimen

Use one of the thin slide holders listed below that cause less interference with the objectives\*.

• U-HRD-4 • U-HLD-4 • U-HLS-4

\* The applicable objectives are ones with 40X or less magnification (except Apo series).

## Using an Oil Immersion Condenser

Use the optional U-SVRO (right-hand knob) or USVLO (left-hand knob) slotted stage which prevents close contact between the specimen and stage.

# Using the Auxiliary Slide Holder

# Applicable revolving nosepieces: U-D7RE, U-D7RES, U-P6RE, U-D6BDRE, U-P5BDRE

When a biological slide glass specimen is placed on the U-HLDT4/U-HRDT4 two-slide holder (thick type), the slide holder may interfere with the objective when the following conditions are combined;

- The objective in use has a short WD (Working Distance);
- The sspecimen being observed is the specimen placed on the rear side of the side holder, and;
- The stage is in a rotated position.

If the slide holder contacts the objective, set the provided auxiliary slide holder ① as shown in the figure and observe only the specimen on the front side of the slide holder.

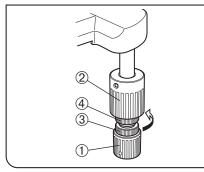


Fig. 27

# 3 Adjusting the X- and Y-Axis Knob Tension (Fig. 27)

- 1. Hold the X-axis knob ① and slide up the Y-axis knob ② up to expose the adjustment knobs.
- 2. Turning the X-axis adjustment knob ③ or Y-axis adjustment knob ④ clockwise (in the direction of the arrow) increases the tension and counterclockwise decreases it.



- If the tension is adjusted to tight, a creaking sound may be heard during stage travel, and the stage stopping accuracy may be imperiled.
- After long hours of use, the stage guide may be deviated and the stage travel range may be decreased.
  - However, this is not malfunction and can be corrected easily as described below.

#### [Treatment]

Horizontal direction: Hold the specimen holder and move the stage guide to the left and right so that it hits the stoppers.

Vertical direction: Hold the upper stage and move it to the front and rear so that it hits the stoppers.

## Stage Feed Knob Rubber Caps (Optional)

When the X- and Y-axis knobs are fitted with the rubber caps, the knobs can be adjusted without slipping and fine adjustment is possible by holding the knobs with a very light force. The knob rubber caps also reduce fatigue after long hours of operation.

The U-SHGT thick type (thickness 5 mm) and U-SHG thin type (thickness 2 mm) knob rubbers are available.

## To attach the knob rubbers:

First fit the larger knob rubber to the Y-axis (upper) knob from below it, then fit the smaller knob rubber to the X-axis (lower) knob from below it.

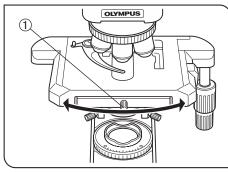


Fig. 28

## 4 Rotating the Stage

- 1. Slightly loosen the stage clamping screw ①.
- 2. The stage can be rotated both clockwise and counterclockwise by the stage clamping screw.

CAUTION A click may be heard and felt during rotation. However, this is due to the construction of the substage and does not indicate a malfunction.

The angle of rotation varies depending on the positions of the X- and Y-axis knobs.

	Angle of Rotation		
	Clockwise	Counterclockwise	
Right hand knobs	230°	20°	
Left hand knobs	20°	230°	


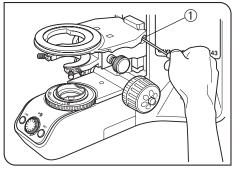


Fig. 29

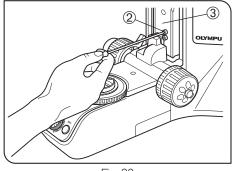

#### Fig. 30

## 5 Adjusting the Stage Height

(Figs. 29 & 30)

- O By lowering the position of the substage, the microscope will accommodate specimens with maximum height of 35 mm. This is useful when observing metallurgical specimens and other thick objects.
- 1. Lower the stage to the lower limit, then remove the stage from the microscope.
- 2. Using the Allen screwdriver, loosen the substage bracket clamping screw ① and remove the substage.
- 3. Turn the coarse adjustment knob and raise the focusing block ③ to where the stopper screw ② on the arm becomes visible.
- 4. Using the Allen screwdriver, loosen and remove the upper stopper screw ②.
- 5. Reattach substage bracket and stage.
- Store the removed stopper screw ② in a safe place so that it will not be lost, if needed again.

# 5-4 Observation Tube

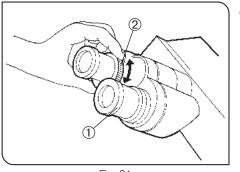


Fig. 31

# Adjusting the Diopter

1

When diopter adjustment rings are provided on both the left and right, it is possible to adjust the optimum vision for both the left and right eyes as described below. The diopter adjustment rings may be provided either on the eyepieces or the observation tube, and there is no problem even when the rings are not provided on both of them.

The diopter adjustment ensures parfocality from low to high magnifications and makes it possible to avoid defocusing even when the objective is switched.

- 1. Set the diopter adjustment rings on both sides to scale "0".
- 2. Engage a high-power objective (40X or so) in the light path, look into the right eyepiece with your right eye, and rotate the coarse and fine adjustment knobs to bring the specimen into focus.
- O not use an oil-immersion objective.
- 3. Engage a low-power objective (10X or so) in the light path, rotate the right diopter adjustment ring ① to bring the specimen into focus. Do not touch the coarse and fine adjustment knobs during the focusing.
- 4. Looking into the left eyepiece with your left eye, rotate the left diopter adjustment ring <sup>(2)</sup> to bring the specimen in focus.
- The above procedure adjusts the diopter with reference to the right eye, but it is also possible to adjust with reference to the left eye. In this case, read the above procedure by inverting "right" and "left."
- If an eyepiece equipped with a diopter adjustment ring is inserted into the side equipped with another diopter adjustment ring of the observation tube, perform the above procedure without moving the diopter adjustment ring of the observation tube from the "0" position.

Fig. 32

# 2 Using the Eye Shades

(Fig. 32)

## When Wearing Eyeglasses

Use with the eye shades in the normal, folded-down position. This will prevent the eyeglasses from being scratched.

## When Not Wearing Eyeglasses

Extend the folded eye shades in the direction of the arrow to prevent extraneous light from entering between the eyepieces and eyes.

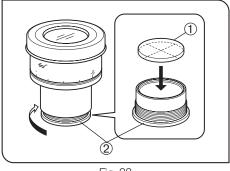


Fig. 33

# **3** Using the Eyepiece Micrometer Disk

(Fig. 33)

When the WHN10X-H (WHN10X) eyepieces are used, an eyepiece micrometer disk can be inserted in one of them. When the eyepiece does not have a diopter adjustment mechanism, however, it is hard to focus on the micrometer disk if the operator has poor eyesight. Should that be the case, adjust the focus with eyeglasses on.

Use 24 mm dia. x 1.5 mm thick micrometer disks.

Following Fig. 33, turn the built-in micrometer-mounting frame @ counterclockwise to remove it from the eyepiece and place a micrometer disk ①.

Screw the micrometer mounting frame back into the eyepiece sleeve. @Be careful not to touch the lens surface with your finger.

The micrometer-mounting frame may be tight for certain micrometer disks. In this case, turn the frame slowly but with a stronger force than usual.

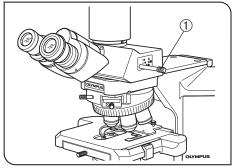


Fig. 34

# 4 Selecting the Light Path of the Trinocular Tube

Slide the light path selector knob to select the desired light path.


Tripper Joy tube	Light path selector knob position			
Trinocular tube	Pushed in	Middle position	Pulled out	
U-TR30-2		20% for binocu-		
U-SWTR-3	100% for bino-	lar eyepieces, 80% for camera		
U-TR30NIR*	cular eyepieces	50% for binocu- lar eyepieces, 50% for camera	100% for camera	
U-TTR2	50% for binocu- lar eyepieces, 50% for camera	100% for binocu- lar eyepieces		

\* With the infrared trinocular tube, infrared observation up to 1000 nm is possible. For details, consult your Olympus representative.

The light path selector knob is located on the right side of the observation tube, but it can be removed and attached on the left side when the U-TTR is used.

After changing the knob position, be sure to tighten the knob firmly with a coin.

(Figs. 35 to 38)

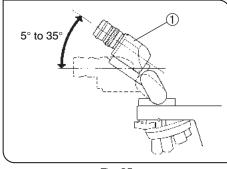



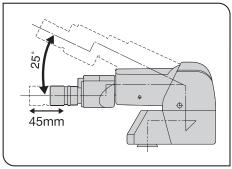


Fig. 36

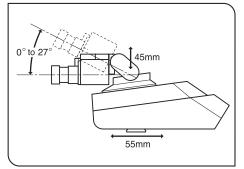


Fig. 37

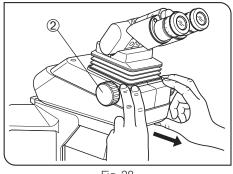


Fig. 38

# 5 Adjusting the Tilt

## With the U-TBI-3 (Fig. 35)

OAdjust the height and tilt of the observation tube to obtain the most comfortable viewing position.

U-TBI-3	5° to 35°
U-TBI-3-CLT	5° to 35°
U-ETBI	0° to 25°
U-TTBI	0° to 25°
U-TTR2	5° to 35°
U-TTLBI	0° to 27°

Holding the binocular section  $\ensuremath{\textcircled{}}$  with both hands, raise or lower it to the desired position.

# CAUTION Never attempt to force the binocular section past the upper or lower stop position. Applying excessive force could destroy the limiting mechanism.

- The U-TBI-3/U-TBI-3-CLT can be used in combination with only one intermediate attachment.
- ◎ If you need photography using the U-TBI-3/U-TBI-3-CLT, use the U-TRU trinocular tube.

## With the U-ETBI/U-TTBI (Fig. 36)

The U-ETBI and U-TTBI are ergonomic observation tubes with normal field, capable of the adjustments of the positioning angle (0° to 25°) as well as front-rear positioning (by 45 mm) of the eyepieces. The U-ETBI is the erect image model and the U-TTBI is the inverted image model, and both models are of the same size.

CAUTION The intermediate attachments that can be combined with the U-TTBI are limited. For details, please contact Olympus.

## With the U-TTLBI (Fig. 37)

The U-TTLBI is a Tilting, Telescopic, Lifting binocular tube with the following adjustment capabilities.

- Eyepiece angle adjustment (0° to 27°)
- Eyepiece up-down position adjustment (45 mm)
- Observation tube front-rear position adjustment (55 mm)

♥When adjusting the front-rear position of the observation tube, hold the base of the observation tube as shown in Fig. 38.

When storing the microscope, increase the angle of the eyepieces so as to prevent the eyepieces from dropping.

◎To adjust the up-down position of the eyepiece, turn the dial ②.

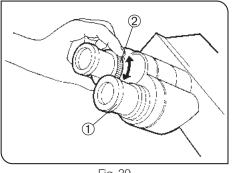


Fig. 39

# 6 Using Eyepieces Incorporating a Micrometer

- When the eyepieces in use incorporate a micrometer, the accuracy of the left-right focusing adjustment (diopter adjustment) can be improved further.
- 1. Looking into the right eyepiece with your right eye, turn the top of the eyepiece ① so that the micrometer in the field of view looks sharpest (Fig. 39).
- 2. Looking into the right eyepiece, rotate the coarse and fine adjustment knobs to bring the specimen and micrometer (cross lines, etc.) in focus.
- 3. Looking into the left eyepiece with your left eye, rotate the diopter adjustment ring <sup>(2)</sup> to bring the specimen into focus.

# 5-5 Condenser

For how to adjust the centering of the condenser, see page 12.

# Compatibility of Objectives and Condensers

	Condenser					
Objective Magnification	Low magnification U-LC	Abbe U-AC2	Achromat/aplanat U-AAC	Swing-out U-SC3	Ultralow magnification U-ULC-2	
1.25X				Applicable (FN 22)		
2X				Applicable (FN 26.5)	Applicable	
4X	Applicable (FN 22)**	Applicable (FN 22)		by moving top lens out of the light path* (FN 26.5)	(FN 26.5)	
10-60X		Applicable	Applicable	Engage top lens in		
100X		(FN 26.5)	(FN 26.5)	light path (FN 26.5)		

\* When using the U-SC3 swing-out condenser together with a 1.25X to 4X objective, fully open the condenser aperture iris diaphragm and use the field iris diaphragm in the base as aperture diaphragm. With the 1.25X to 2X objectives, the surroundings of the field of view may become dark.

To obtain better illumination in camera recording using a 1.25X to 4X objective, use of the U-ULC-2 is recommended.
\*\*When the U-LC is used with an objective with an NA of 1.0 or more, the NA of the lighting becomes insufficient.
When the U-LC is used, it is required to attach the auxiliary lens (page 34).

# 5-6 Immersion Objectives

CAUTION B

Be sure to use the provided Olympus Immersion oil.

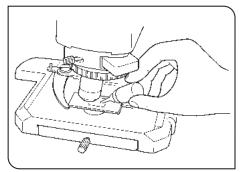


Fig. 40

# Using Immersion Objectives

(Fig. 40)

- 1. Focus on the specimen with objectives in the order of lower-power to higher-power ones.
- 2. Before engaging the immersion objective, place a drop of provided immersion oil onto the specimen at the area to be observed.
- 3. Turn the revolving nosepiece to engage the immersion objective, then focus using the fine adjustment knob.



Since air bubbles in the oil will affect the image quality, make sure that the oil is free of bubbles.

- a. To check for bubbles, remove the eyepiece and fully open the field and aperture iris diaphragms, then look at the exit pupil of the objective inside the observation tube. (The pupil should appear round and bright.)
- b. To remove bubbles, turn the revolving nosepiece to repeatedly defocus and refocus the oil immersion objective.
- If the condenser engraving shows a numerical aperture (NA) of 1.0 or higher, the number applies only when oil is applied between the slide glass and the top surface of the condenser. When oil is not present, the NA is about 0.9.
- 4. After use, remove immersion oil from the objective front lens by wiping with gauze slightly moistened with absolute alcohol.

#### CAUTION


#### Caution in use of immersion oil

If immersion oil enters your eyes or contacts with your skin, immediately take the following treatment.

- Eyes: Rinse with fresh water (for 15 minutes or more).
- Skin: Rinse with water and soap.

If the appearance of the eyes or skin is altered or pain persists, immediately see your doctor.

# 5-7 Objectives with Correction Collar

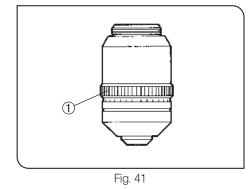

If the thickness scale setting of the objective with collection collar does not match the cover glass thickness, the objective cannot manifest the full performance. To prevent this, be sure to perform the following adjustment when using an objective with collection collar.

#### Adjustment Procedure

- If the cover glass thickness is known, set the correction collar ① to that scale value. (Fig. 41)
- If the cover glass thickness is unknown, adjust the correction collar ① and fine adjustment knob alternately until the positioning with the highest contrast is obtained.
- Be careful not to touch the correction collar ① when turning the revolving nosepiece.


© Use a trinocular observation tube such as the U-TR30-2 or U-SWTR-2 when shooting the video or digital camera picture of the microscope images.

The trinocular tube accepts a camera adapter (certain TV adapters necessitate a camera mount adapter).

Be sure to adjust the parfocality before using a camera adapter. Otherwise, the focusing of the camera image will not
 match that of the image observed through eyepieces. For the parfocality adjustment method, refer to the instruction
 manual for the camera adapter in use.

For details, refer to the instruction manuals for these accessories.


## System Chart

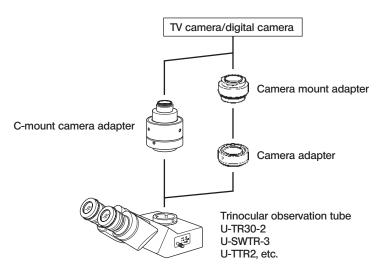

# 2 Selecting the Camera Adapter Magnification

The image pickup area is determined by the size of the image sensor used in the TV camera or digital camera and the magnification of the camera adapter.

The following illustrations show the image pickup areas of camera adapters when the SWH10X-H eyepieces with a FN of 26.5 are used (the FN of the WHN10X eyepieces is 22).

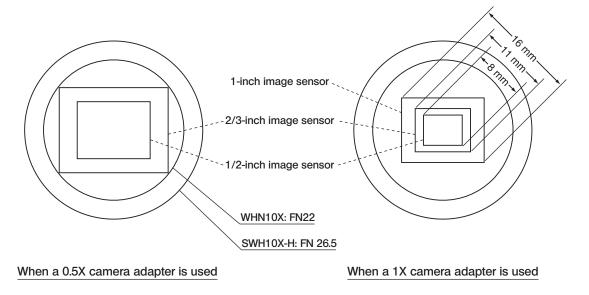

# TROUBLESHOOTING GUIDE

7

Under certain conditions, performance of the unit may be adversely affected by factors other than defects. If problems occur, please review the following list and take remedial action as needed. If you cannot solve the problem after checking the entire list, please contact your local Olympus representative for assistance.

Problem	Lamp	Cause	Remedy	Page
Bulb does not light.	LED	Power is not ON.	Set the main switch to " I " (ON).	10
	Common	Power cord is unplugged.	Plug power cord into the power outlet.	37
	Common	Connectors are connected improperly.	Connect connectors securely.	37
	Halogen	Bulb is not attached.	Attach a halogen bulb.	34
	Halogen	Bulb is burned out.	Replace bulb.	34
	Halogen	The TL4 is not ON.	Set the main switch of the TL4 to "I" (ON).	37
Bulb is blown immediately.	Halogen	The bulb is not a designated one.	Replace bulb with a designated one.	1
Field of view is dark.	LED	The LIM setting is inappropriate.	Set LIM again.	13
	Common	Aperture and field iris diaphragms are not opened wide enough.	Adjust them to proper sizes.	12
	Common	Light path selector knob is set to the rôp position.	Set the knob to the ôd or a position.	22
	Common	Condenser is not properly centered.	Center the condenser.	12
	Common	Condenser height is too low.	Adjust correctly.	24
	Halogen	ND filter is engaged.	Disengage the ND filters.	15
Brightness cannot be varied.	LED	LIM is not ON.	Switch the LIM function OFF.	13
	Halogen	LED brightness adjustment knob is rotated.	Use the brightness adjustment knob of the TL4.	37
	Common	Condenser is not properly centered.	Center the condenser.	12
	Common	Condenser height is too low.	Adjust correctly.	24
Field of view is obscured or not evenly illuminated.	Common	Light path selector knob is set in an intermediate position.	Set to the desired click position.	22
	Common	Objective is not correctly engaged/	Make sure that the revolving nosepiece clicks properly into place.	12
	Common	Revolving nosepiece is not attached properly.	Push the slide dovetail all the way until it is stopped.	35
	Common	Condenser is not attached properly.	Re-attach it .	36
	Common	An objective that falls outside of the condenser's illumination range is used.	Use a condenser to match the purpose.	24
	Common	Condenser is not properly centered.	Center the condenser.	12
	Common	Field iris diaphragm is stopped down too far.	Open the field iris diaphragm until it circumscribes the field.	12
	Halogen	Bulb is not mounted correctly.	Push the pins of halogen bulb all the way until the stop position.	34

Problem	Lamp	Cause	Remedy	Page
Dirt or dust is visible in the field	Common	Dirt/dust on the eyepieces.	Clean thoroughly.	
of view.	Common	Dirt or dust on condenser top lens.		3
	Common	Dirt/dust on the specimen.		
Visibility is poor • Image is not poor.	Common	A non-UIS/UIS2 objective is used.	Use only UIS2/UIS series objectives with this microscope.	32
Contrast is poor.	Common	Condenser height is too low.	Adjust correctly.	24
<ul><li>Details are indistinct.</li><li>Image glares.</li></ul>	Common	Aperture iris diaphragm is not opened enough.	Open aperture iris diaphragm.	12
	Common	Revolving nosepiece is not at- tached properly.	Push the slide dovetail all the way until it is stopped.	35
	Common	Correction collar on correction collar-equipped objective is not properly adjusted.	While focusing, turn the correction collar to find the best position.	25
	Common	Front lens of objective is dirty.	Clean objective.	3
	Common	Immersion oil is not being used with an oil immersion objective.	Use immersion oil.	25
	Common	Immersion oil contains bubbles.	Remove the bubbles.	25
	Common	Recommended immersion oil is not used.	Use an Olympus-designated immersion oil.	25
	Common	Dirt/dust on specimen.	Clean it.	3
	Common	Dirt/dust on condenser top lens.		
One side of image is blurred.	Common	Objective is not correctly engaged in light path.	Make sure that revolving nosepiece clicks into place correctly.	12
	Common	Stage is not correctly mounted.	Re-attach it.	18
	Common	Revolving nosepiece is not at- tached properly.	Push the slide dovetail all the way until it is stopped.	35
	Common	Specimen is mounted upside down.	Mount specimen correctly.	18
Image appears to waver.	Common	Revolving nosepiece is not attached properly.	Push the slide dovetail all the way until it is stopped.	35
	Common	Objective is not correctly engaged in light path.	Make sure that revolving nosepie- ce clicks into place correctly.	12
	Common	Condenser is not properly centered.	Center the condenser.	12
Coarse adjustment knob is hard to turn.	Common	Tension adjustment ring is tightened excessively.	Loose ring.	16
	Common	You are trying to raise stage without coarse adjustment knob while pre-focusing lever is kept locked.	Unlock pre-focusing lever.	17
<ul><li>Image cannot be focused:</li><li>Objective drifts down by itself.</li><li>Fine adjustment knob slips.</li></ul>	Common	Tension adjustment ring is too loose.	Tighten ring.	16

Problem	Lamp	Cause	Remedy	Page
Coarse adjustment will not go all the way up.	Common	Pre-focusing lever is locked at a low position.	Unlock pre-focusing lever.	17
Coarse adjustment will not go all the way down.	Common	Condenser holder is too low.	Raise condenser holder.	36
Objective makes contact with specimen before focus is obtained.	Common	Specimen is mounted upside down.	Mount specimen correctly.	18
Field of view of one eye does not	Common	Interpupillary distance is incorrect.	Adjust interpupillary distance.	11
match that of the other.	Common	Diopter adjustment is incorrect.	Adjust diopter.	21
	Common	Different eyepieces are used on left and right.	Change one eyepiece to match the other so that both sides are the same type.	32
	Common	Your view is not accustomed to microscope observation.	Upon looking into eyepieces, try looking at overall field before concentrating on specimen ran- ge. You may also find it helpful to look up and into distance for a moment before looking back into microscope.	21
Image shifts when you touch stage.	Common	Stage is not properly mounted.	Clamp stage.	18
X- and Y-axis knobs are too tight or too loose.	Common	Tension of X- and Y-axis knobs is too high or too low.	Adjust tension.	19
Stroke has reduced.	Common	Stage guide is deviated.	Correct deviation as described.	19



Item	Specification							
1. Optical system	UIS2 (UIS) optical system (featuring infinity correction)							
2. Illumination	Built-in transmitted Koehler illumination         FN (Field Number): 26.5 (super-widefield compatible)         LED illuminator (continuously variable): U-LHLEDC         Power consumption: 0.5 A         Optional       6V 30W halogen bulb (continuously variable)         Power supply unit TL4: 100-120/220-240 V ∼ , 0.85/0.45 A, 50/60 Hz.         Power consumption: 45 W							
3. Focusing mechanism	Stage up-down movement by roller guide (rack & pinion) Stroke per rotation: 0.1 mm (fine), 17.8 mm (coarse) Full stroke range: 25 mm Upper coarse adjustment limit stopper Tension adjustment on coarse focus adjustment knob.							
4. Revolving nosepiece	Туре	BX43	-5RES	U-D	06RE			
		5-position revolvir readout	ng nosepiece with function	Universal 6-position	revolving nosepiece			
	Attachable module			ght DIC slider light analyzer				
5. Observation tube	Туре	U-BI30-2	U-TBI-3/U-TBI3-CLI	U-TR30-2	U-SWTR-3			
		Widefield binocular	Widefield, tilting binocular	Widefield, trinocular	Super- widefield, trinocular			
	Field No.		22		26.5			
	Tube inclination	30° 5-35° (continuously variable)		30°				
	Interpupillary distance adjustment	50-76 mm						
	Light path selection	Not av	ailable	3 steps: ① Binocular 100% ② Bi 100%, Cam 80% ③ Camera 100%				
6. BX stage	Туре	U-SVRB-4		U-SVLB-4				
		Coaxial knobs	on bottom right.	Coaxial knobs on bottom left.				
		Rectangular ceramic-coated, wire-driven stage						
	Size	156 mm (D) x 191 mm (W)						
	Movement mechanism	X- and Y-axis knob with adjustable tension. Movement range: 52 mm in vertical (Y) direction, 76 mm in horizontal (X) direction.						
	Slide holder (single-slide holder)	U-HLST-4 (Thick type) U-HLS-4 (Thin type)						
	Slide holder (double-slide holder)		pening, thick type) pening, thin type)	U-HRDT-4 (Right-opening, thick type) U-HRD-4(Right-opening, thin type)				

Item		Specification							
7. Condenser	Туре	U-LC	U-AC2	U-SC3	U-AAC				
Max. N.		Low magnification	Abbe	Swing-out	Achromat/ Aplanat				
		0.75	1.10	0.9	1.40				
	Applicable objective powers	2X - 60X (Widefield: FN 22)	4X (Widefield: FN 22) 10X - 100X (Super widefield: FN 26.5)	1.25X (Widefield: FN 22) 2X - 100X (Super widefield: FN 26.5)	10X - 100X (Super widefield: FN 26.5)				
8. Operating environment	<ul> <li>Ambient ter</li> <li>Maximum r 70% at 34°</li> <li>Supply volta</li> <li>Pollution de</li> </ul>	<ul> <li>Indoor use.</li> <li>Altitude: Max. 2000 meters</li> <li>Ambient temperature: 5° to 40°C (41° to 104°F)</li> <li>Maximum relative humidity: 80% for temperatures up to 31°C (88°F), decreasing linearly through 70% at 34°C (93°F), 60% at 37°C (99°F), to 50% relative humidity at 40°C (104°F).</li> <li>Supply voltage fluctuations: ±10%</li> <li>Pollution degree: 2 (in accordance with IEC60664-1)</li> <li>Installation/Overvoltage category: II (in accordance with IEC60664-1)</li> </ul>							
9. Rating		Input rating: 5 V, 2.5 A AC adapter SA115B-05V: 100V-240Vac 50/60 Hz, 0.4 A							

**OPTICAL CHARACTERISTICS** (UIS2 Series)


#### - UIS series objectives not listed here can also be combined with this microscope. -

The following table shows the optical characteristics of combinations of eyepieces and objectives. The figure on the right shows the performance data engraved on the objectives.

#### (NOTE )

Y

Refer to the latest catalogue or consult your local Olympus representative for the updated information on the eyepieces and objectives that can be combined with this microscope.

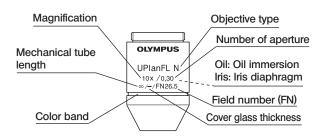

Optical character		Magnifi-		W.D.		Reso-	Eyepiece WHN10X (FN22)				
Objective Series	Marking	cation	N.A.	(mm)	thick- ness (mm)	lution (µm)	Total mag.	Depth of focus (µm)	Field of view	Remark	
PLN (PH) Plan Achromat (FN 22) PH: Phase contrast	PlanN (Ph)	2X 4X 10X (Ph) 20X (Ph) 40X (Ph) 50XOI 100XO (Ph)	0.06 0.1 0.25 0.4 0.65 0.5-0.9 1.25	5.8 18.5 10.6 1.2 0.6 0.2 0.15	- 0.17 0.17 0.17 -	5.59 3.36 1.34 0.84 0.52 0.37 0.27	20X 40X 100X 200X 400X 500X 1000X	560.1 175.0 28.0 9.27 3.04 1.7 0.69	11.0 5.5 2.2 1.1 0.55 0.44 0.22	Oil immersion/Iris Oil immersion	
UPLFLN (PH) Plan Semi Apochromat (FN 26.5) PH: Phase contrast	UPlanFLN (Ph)	4X (Ph) 10X2 (Ph) 20X (Ph) 40X (Ph) 40XO 60X 60XOI (Ph) 100XO2 (Ph) 100XOI2	0.13 0.3 0.5 0.75 1.3 0.9 0.65-1.25 1.30 0.6-1.30	17.0 10.0 2.1 0.51 0.2 0.2 0.12 0.2 0.2 0.2	- 0.17 0.17 0.17 0.17 0.17 0.17 0.17	2.58 1.12 0.67 0.45 0.26 0.37 0.27 0.26 0.26	40X 100X 200X 400X 400X 600X 600X 1000X 1000X	127.2 22.4 7.0 2.52 1.27 1.5 0.98 0.66 0.66	5.5 2.2 1.1 0.55 0.37 0.37 0.22 0.22	Oil immersion Correction collar Oil immersion/Iris Oil immersion Oil immersion/Iris	
UPLSAPO Plan Apochromat (FN 26.5)	UPlanSApo	4X 10X2 20X 20XO 40X2 60XW 60XO 100XO	0.16 0.4 0.75 0.85 0.95 1.2 1.35 1.4	13.0 3.1 0.6 0.17 0.18 0.28 0.15 0.13	- 0.17 0.17 - 0.17 0.17 0.17 0.17	2.10 0.84 0.45 0.39 0.35 0.28 0.25 0.24	40X 100X 200X 200X 400X 600X 600X 1000X	99.6 15.9 4.29 3.5 1.9 1.03 0.89 0.59	5.5 2.2 1.1 1.1 0.55 0.37 0.37 0.22	Oil immersion Correction collar Water immersion Oil immersion Oil immersion	
PLAPON Plan Apochromat (FN 26.5)	PlanApoN	1.25X 2X 60XO	0.04 0.08 1.42	5.0 6.2 0.15	- - 0.17	8.39 4.19 0.24	12.5X 20X 600X	1326.8 398.3 0.83	17.6 11.0 0.37	Oil immersion	
UApoN Apochromat (FN 22)	UApoN	20XW340 40XW340 40XO340	0.7 1.15 1.35	0.35 0.25 0.1	0.17 0.13-0.25 0.17	0.48 0.29 0.25	200X 400X 400X	4.08 1.29 1.2	1.1 0.55 0.55	Water immersion Water immersion/ Correction collar Oil immersion	




### 10-1 Assembly Diagram

The diagram below shows the sequence of assembly of the various modules. The numbers indicate the order of assembly.


Assembly steps enclosed in 
will be detailed on the subsequent pages.

The module numbers shown in the following diagram are merely the typical examples. For the modules with which the module numbers are not given, please consult your Olympus representative or the catalogues.

## CAUTION When assembling the microscope, make sure that all parts are free of dust and dirt, and avoid scratching any parts or touching glass surfaces.

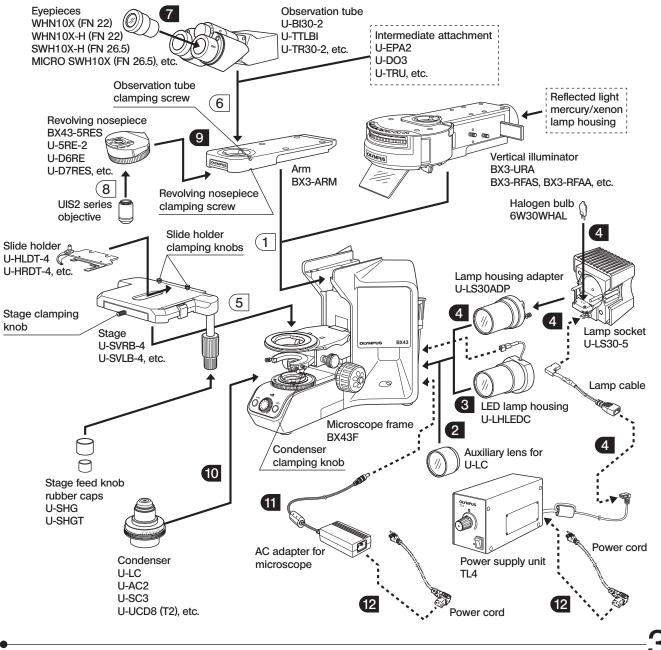

### **10-2 Detailed Assembly Procedures**

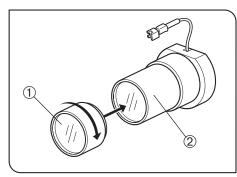

# Attaching the Auxiliary Lens for the U-LC (Fig. 42)


- ♥It is not necessary to attach the auxiliary lens if the condenser in use is other than the U-LC.
- 1. Attach the auxiliary lens ① provided with the condenser into the tip of the LED lamp housing C (U-LHLEDC) ② or lamp housing adapter (U-LS30ADP) by screwing the lens in the direction of the arrow.



2

#### Be careful not to touch the lens with your finger.

Fig. 42

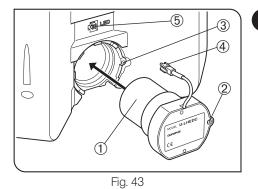

#### 3 Attaching the LED Lamp Housing C

 Fit the LED lamp housing ① into the mount hole on the rear of the microscope by aligning the clamping screw ② and screw hole ③.
 Using the Allen screwdriver, tighten the clamping screw ②.

(Fig. 43)

3. Insert the connector ④ of the lamp housing into the connector ⑤ on the microscope.

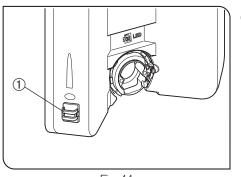


Fig. 44

### 4 Attaching the Halogen Lamp Socket (Fig. 44)

#### Installing the U-LS30ADP LS30 Adapter (Fig. 44)

- 1. Peel off the backing of double-side adhesive tape on the cable holder 1 and attach the holder on the specified position.
- The area (enclosed with broken lines) on the outer side is for use in attaching the cable holder for another option.
- 2. Attach the adapter in the same way as steps 1 and 2 in the LED lamp housing C attaching procedure described above.

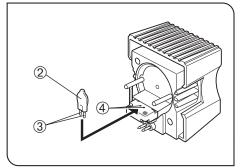


Fig. 45

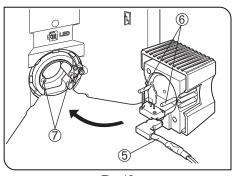


Fig. 46


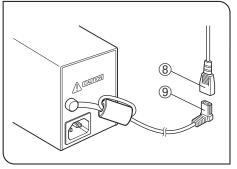


Fig. 47

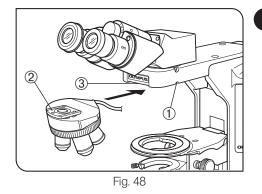


Installing the Halogen Bulb (Fig. 45)

#### Use only the designated bulb 6V30WHAL (PHILIPS 5761).

1. Holding the bulb ② a piece of gauze, insert the bulb pins ③ straight and fully into the pin holes ④ on the lamp socket.

CAUTION To

To prevent reduced bulb life or cracking, do not touch the bulb with bare hands. If fingerprints are accidentally left on the bulb, wipe the bulb with a soft cloth.

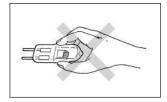

2. Insert the lamp cable (5) in the direction shown in the figure 46.
3. Aligning the guide pins (6) with the guide pin holes (7) at the rear of the microscope frame, gently push the lamp socket all the way into place.



Caution for Bulb Replacement During Use or Right After Use

The bulb and the lamp socket are areas near these will be extremely hot during and right after use.

Set the main switch to "  $\mathbf{O}$  " (OFF), disconnect the power cord from the wall outlet, then allow the old bulb and lamp socket to cool before replacing the bulb with a new of the designated type.


4 Pass the lamp cable through the cable holder (Fig. 44).

#### Connecting the Lamp Cable (Fig. 47)

Connect the cable connector (18) to the connector (19) of the TL4 power supply unit.

#### 8 Attaching the Revolving Nosepiece (Fig. 48)

- 1. Rotate the coarse adjustment knob to lower the stage.
- 2. Loosen the revolving nosepiece clamping screw ① slightly. Do not loosen it too much because doing so makes it impossible to attach the revolving nosepiece.
- Hold the revolving nosepiece so that the mount dovetail <sup>(2)</sup> is horizontal and insert it all the way into the dovetail groove of the microscope arm <sup>(3)</sup>.
- 4. Tighten the revolving nosepiece clamping screw 1.
- When attaching or detaching a revolving nosepiece with a cable, be sure to disconnect the cable from the revolving nosepiece in advance.

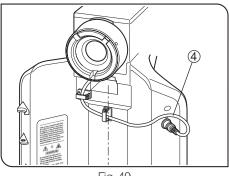

#### Use only coded 5 position nosepiece for BX43 (BX43-5RES)

- Obstribute its cable ④ toward the rear panel by passing it below the arm, and connect the cable to the connector no the rear panel.

Fig. 49

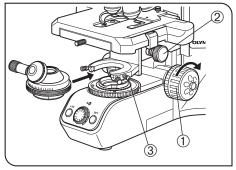


Fig. 50


#### 9 Attaching the Condenser

(Fig. 50)

- Turn the coarse adjustment knob ① to raise the stage to its upper limit.
   Turn the condenser height adjustment knob ② to lower the condenser holder to its lowest position.
- 3. Fully loosen the condenser clamping knob 3.
- 4. Position the condenser so that the aperture scale is located in front. Gently slide the condenser along the dovetail all the way until it is stopped.

<sup>©</sup>When attaching the U-SC3 swing-out condenser, align the positioning pin at the back of the condenser with the groove in the condenser sleeve.

- When attaching a swing-out condenser with a positioning pin at the back such as the U-SC3, attach by aligning the positioning pin with the groove in the condenser sleeve.
  - When attaching a condenser with swing-out top lens, swing out the top lens out of the way before inserting the condenser.
  - 5. Tighten the condenser clamping knob and gently raise the condenser holder to its highest position.

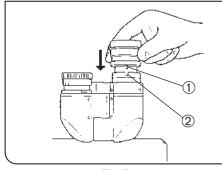



#### 10 Attaching the Eyepieces


(Fig. 51)

Gently insert the eyepieces all the way into the eyepiece sleeves.

- When using the U-BI30-2 binocular tube, eyepieces with a positioning pin cannot be attached because the U-BI30-2 does not have the positioning notch.
  - When using an eyepiece with micrometer, attach it to the righthand eyepiece sleeve.

When doing so, make sure that the eyepiece positioning pin  ${\rm (}0$  fits into the notch  ${\rm (}2$  on the eyepiece sleeve.

• The super-widefield trinocular observation tube is equipped with a positioning notch on both the left and right eyepiece sleeves. Make sure that the positioning pins on both eyepieces fits into the respective notches.

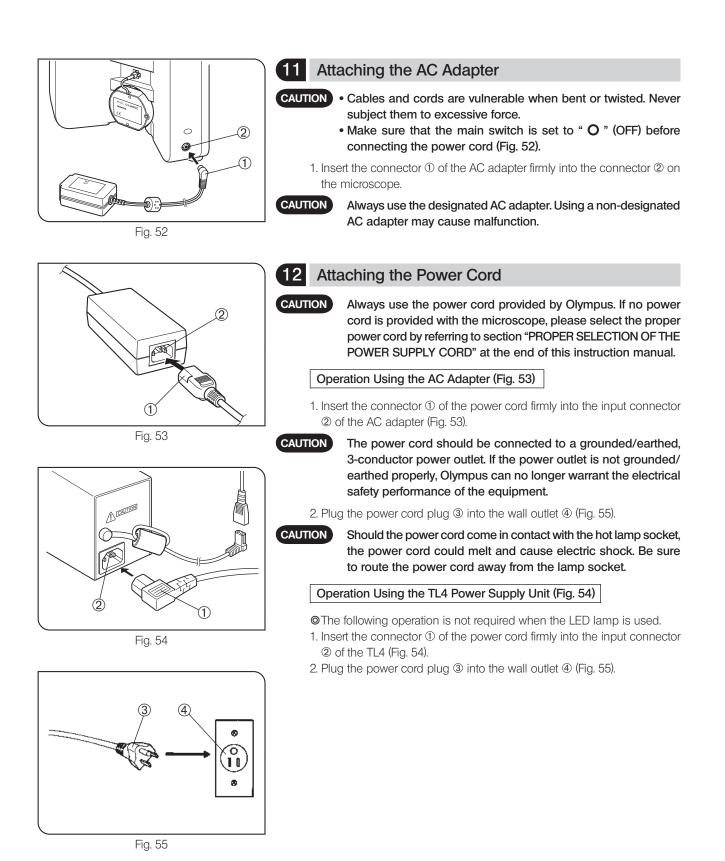

- For safe use of the lamp socket, we recommend performing the following inspection periodically (every time you replace the lamp bulb and at least every 6 months).
- The table below identified the check items to be observed. Put (X) if not applicable or ( $\checkmark$ ) if applicable.
- If there is any ( $\checkmark$ ) mark noted, immediately stop use of the product, and consult Olympus for detailed inspections or replace the lamp socket.
- If you detect an abnormality other than that listed below or with other Olympus product, also stop the use of the product and contact Olympus for detailed inspections.
- •Note that the service, replacement and detailed inspections are charged after expiration of the warranty period.

If you have any questions, please contact Olympus.

	Check results (Date)			
Check items		/	/	/
1. More than 8 years have passed since original purchase or the total power ON time exceeds 20,000 hours.				
2. Lamp does not light sometimes even though the main switch is set to on.				
3. Illumination flickers when you move the lamp cable or lamp socket.				
4. Scorching or burning odor is produced during use.				
5. Illumination still flickers after replacement with a new lamp bulb.				
6. Deformation, backlash, or looseness, etc. when you assemble the lamp socket.				
7. Extreme discoloration of the lamp socket connection terminal or lamp socket lamp bulb mount.				
8. Discoloration, deformation or cracking of the lamp socket.				
9. Melting, crack, deformation or solidification of the lamp cable or a wiring part.				
10. Increased frequency of servicing compared to similar devices put into use at the same time as the lamp socket.				

\* When the Check Result columns become insufficient, copy this sheet.

### ■ PROPER SELECTION OF THE POWER SUPPLY CORD

If no power supply cord is provided, please select the proper power supply cord for the equipment by referring to "Specifications" and "Certified Cord" below:

## CAUTION: In case you use a non-approved power supply cord for Olympus products, Olympus can no longer warrant the electrical safety of the equipment.

#### **Specifications**

Voltage Rating	125V AC (for 100-120V AC area) or, 250V AC (for 220-240V AC area)
Current Rating	6A minimum
Temperature Rating	60°C minimum
Length	3.05 m maximum
Fittings Configuration	Grounding type attachment plug cap. Opposite terminates in molded-on IEC configuration appliance coupling.

#### Table 1 Certified Cord

A power supply cord should be certified by one of the agencies listed in Table 1, or comprised of cordage marked with an agency marking per Table 1 or marked per Table 2. The fittings are to be marked with at least one of agencies listed in Table 1. In case you are unable to buy locally in your country the power supply cord which is approved by one of the agencies mentioned in Table 1, please use replacements approved by any other equivalent and authorized agencies in your country.

Country	Agency	Certification Mark	Country	Agency	Certification Mark	
Argentina	IRAM	æ	Italy	IMQ	$\odot$	
Australia	SAA	$\Delta$	Japan	JET, JQA, TÜV, UL-APEX/MITI	(PS),	
Austria	ÖVE	<b>Øve</b>	Netherlands	KEMA	KEUR	
Belgium	CEBEC	<b></b>	Norway	NEMKO	N	
Canada	CSA	<u>ج</u>	Spain	AEE	Ø	
Denmark	DEMKO	D	Sweden	SEMKO	\$	
Finland	FEI	F	Switzerland	SEV	(† S	
France	UTE	(S)	United Kingdom	ASTA BSI	€, 🕏	
Germany	VDE	ĎÈ	U.S.A.	UL	(L	
Ireland	NSAI	ø		· · · · · · · · · · · · · · · · · · ·		

#### Table 2 HAR Flexible Cord

#### APPROVAL ORGANIZATIONS AND CORDAGE HARMONIZATION MARKING METHODS

Approval Organization	Printed or Embossed Harmoni- zation Marking (May be located on jacket or insulation of internal wiring)		Alternative Marking Utilizing Black-Red-Yellow Thread (Length of color section in mm)			
			Black	Red	Yellow	
Comite Electrotechnique Belge (CEBEC)	CEBEC	<har></har>	10	30	10	
Verband Deutscher Elektrotechniker (VDE) e.V. Prüstelle	<vde></vde>	(HAR)	30	10	10	
Union Technique de l'Electricite' (UTE)	USE	(HAR)	30	10	30	
Instituto Italiano del Marchio di Qualita' (IMQ)	IEMMEQU	(HAR)	10	30	50	
British Approvals Service for Electric Cables (BASEC)	BASEC	<hr/>	10	10	30	
N.V. KEMA	KEMA-KEUR	(HAR)	10	30	30	
SEMKO AB Svenska Elektriska Materielkontrollanstalter	SEMKO	<har></har>	10	10	50	
Österreichischer Verband für Elektrotechnik (ÖVE)	<öve>	<har></har>	30	10	50	
Danmarks Elektriske Materialkontroll (DEMKO)	(DEMKO)	<har></har>	30	10	30	
National Standards Authority of Ireland (NSAI)	(NSAI)	<har></har>	30	30	50	
Norges Elektriske Materiellkontroll (NEMKO)	NEMKO	<har></har>	10	10	70	
Asociacion Electrotecnica Y Electronica Espanola (AEE)	(UNED)	<har></har>	30	10	70	
Hellenic Organization for Standardization (ELOT)	ELOT	(HAR)	30	30	70	
Instituto Portages da Qualidade (IPQ)	np	(HAR)	10	10	90	
Schweizerischer Elektro Technischer Verein (SEV)	SEV	(HAR)	10	30	90	
Elektriska Inspektoratet	SETI	(HAR)	10	30	90	

Underwriters Laboratories Inc. (UL) Canadian Standards Association (CSA) SV, SVT, SJ or SJT, 3 X 18AWG

SV, SVT, SJ or SJT, 3 X 18AWG

# MEMO

# MEMO





OLYMPUS CORPORATION Shinjuku Monolith, 3-1, Nishi Shinjuku 2-chome, Shinjuku-ku, Tokyo, Japan OLYMPUS EUROPA HOLDING GMBH Wendenstrasse 14-18, 20097 Hamburg, Germany OLYMPUS AMERICA INC. 3500 Corporate Parkway, Center Valley, Pennsylvania 18034-0610, U.S.A. OLYMPUS SINGAPORE PTE LTD. 491B River Valley Road, #12-01/04 Valley Point Office Tower, Singapore 248373 OLYMPUS AUSTRALIA PTY. LTD. 31 Gilby Road, Mount Waverley, VIC., 3149, Melbourne, Australia OLYMPUS LATIN AMERICA, INC. 5301 Blue Lagoon Drive, Suite 290 Miami, FL 33126, U.S.A.

#### Distributor

**Tengrant Inc.** Building 1,704/1199 No. Xing Mei Road,Shanghai,China Tel.: +86(021)5428 6005 Fax: +86(021)5428 6005-8002 www.tengrant.com a88888@sina.cn